

---

## Using LabQuest 2 with an iPad

---

### Switch on LabQuest 2 (it takes a while)

Plug in probe to Channel 1

→ Home page

→ Connections

WiFi ON

Name: LabQuest

Data sharing: ON

Deselect Allow connected devices to start/stop collection

(Otherwise students could switch off the collection by the LabQuest)

→ OK

Network Settings symbol (top right)

Scroll down the networks

→ Create a network

Give it a name, use the number on device (e.g. Labquest 5)

Create

OK

QR Code displays in the box

### Switch on iPad

Open settings

Switch Wifi ON

In Wifi

Chose a network

Select LabQuest...

(WiFi should now say LabQuest... next to it)

### Open Graphical Analysis

→ Create an experiment

Data sharing

→ Specify source

Click **Camera**

→ Scan QR Code

(make sure the camera is enabled on the iPad, it should ask you but if not it is in settings).

Click **Connect**

Connection made with LabQuest...

### Back to Labquest

→ Home

→ LabQuest App

Set Mode/ Frequency of measurement

Press record tab (green arrow)

Graph appears in real time on the iPad

You can add a trend line, curve fit etc.

### Edit

→ Waste bin

→ Delete experiment

Terminate network in Network Wifi connection preferences

Graphical Analysis also exists for Smart phones.